

## 1. Description

Seplife® 6AG IDA/90 is a metal affinity chromatography resin based on 6% highly cross-linked agarose functionalized with iminodiacetic acid (IDA).

- Seplife® 6AG IDA/90 can be loaded with a range of divalent metals such as Ni, Co, Cu, Fe, Zn showing high affinity for His-tagged proteins.
- Suitable for small scale and large scale purifications of His-tagged proteins from other proteins in the feed.
- Seplife® 6AG IDA/90 resin provides high dynamic binding capacity toward His-tag proteins upon loading with divalent metal.
- High stability to CIP (cleaning in place) up to 1.0M NaOH.
- Regulatory Support File (RSF) documentation is available for Seplife® 6AG IDA/90.

Seplife® 6AG IDA/90 is based on highly cross-linked (6%) agarose functionalized with IDA and ready to be loaded with a divalent metal of choice for affinity binding to His-tag proteins and has a particle size range of 45-165 micron.

## 2. Properties

Product	Seplife® 6AG IDA/90
Appearance	White spherical beads
Type	Metal affinity chromatography resin (IDA)
Matrix	6% cross-linked agarose
Ligand	Iminodiacetic acid
Particle size range ( μ m)	45-165
pH stability	3-13 (operational), 2-14 (CIP)
Chemical Stability	Stable in all common aqueous buffers; 1M sodium hydroxide; 8M urea; 6M guanidine hydrochloride; 70% ethanol.
Flow rate* (cm/h)	Max 750cm/h
Shipped as	20% ethanol slurry

\*Testing conditions: Chromatography column 16mm×200mm; column bed height 20cm; temperature 25° C; mobile phase water.

### 3. Instructions

Pretreatment: metals can be chelated according to actual needs. After the metal is chelated, use the media according to the following procedures.

#### 3.1 Column packing

Column packing should be done according to standard operating procedures. It is important to ensure that each material is at its working temperature, and when possible, the chromatography media may be degassed before column packing.

#### 3.2 Equilibration

Equilibrate the column with an appropriate 2-5 column volume buffer. Ensure the conductivity and pH of the effluent are exactly the same as the buffer.

#### 3.3 Sample feeding

1. The sample is generally dissolved in the equilibration buffer solution at pH 6-8, and increasing the pH of the loading buffer may increase the protein loading capacity.
2. The buffer should not contain metal chelating agents such as EDTA and citrate, and it is best to avoid reducing agents such as mercaptoethanol and DTT.
3. Commonly used buffers include 10-100 mmol/L sodium phosphate, 20-200 mmol/L Tris-HCl, etc.
4. Generally, 0.15-0.5mol/L NaCl should be added to the buffer to eliminate ion exchange interactions.
5. When using the metal chelating agarose media for the first time, it is recommended to use 50mmol/L PBS (50 mmol/L NaH<sub>2</sub>PO<sub>4</sub>, 0.5 mol/L NaCl, pH 7.4) as the initial buffer.

#### 3.4 Elution

Typically, the following methods can be used for protein elution:

1. Lower pH: most proteins will be eluted at pH 6-4 (also at pH 3-4), the buffer can be sodium acetate, and phosphate buffer systems.
2. Competitive elution: linearly increase or one-step increase the concentration of competing substances with affinity to metal ions (such as 0-1.0 mol/L imidazole, 0-50 mmol/L histidine, 0-2 mol/L NH<sub>4</sub>Cl).
3. Chelating agents: Chelating agents such as EDTA and EGTA can bind to metal ions and release the proteins. However, this method cannot separate different proteins and will remove the metal from the resin.

#### Note:

1. When using for the first time, if the concentration required for elution is not certain, it is recommended to add 10mmol/L, 20mmol/L, 50mmol/L, 100mmol/L, 200mmol/L, 500mmol/L or more imidazole to the initial

buffer solution, from low to high concentration, to elute and collect the recombinant proteins respectively, and then identify the elution results by SDS-PAGE electrophoresis and other methods. If conditions permit, linear imidazole gradient elution can be performed to determine better elution conditions.

2. Imidazole is alkaline, and the pH needs to be adjusted using HCl after the corresponding buffer solution is prepared.

3. Elution by lowering pH and with chelating agent will cause the metal ions to be removed from the resin. The metal ions will need to be re-chelated to the resin before the next use.

For all the above elution methods, 150-500mmol/L NaCl should be added to the buffer to eliminate any ion exchange interactions.

### 3.5 Regeneration

1. After multiple uses or when the chelated metal ions need to be replaced, the metal must be stripped of the chromatographic media and the resin regenerated.

Metal removal method: First rinse the column with 5-10 CV of distilled water, then rinse the column with 5-10 CV of 100mmol/L EDTA, and finally use 2-3 CV of 0.5mol/L NaCl to wash off residual EDTA.

2. Columns that have been used many times generally require cleaning after metal removal.

Cleaning method: reverse wash the column with 0.1-1.0 mol/L NaOH, and keep it at 50 cm/h for 1-2 hours, which should remove the strongly bound impurities.

3. Metal ions need to be re-chelated after washing.

Chelating method: First, use 2-5 CV of distilled water to fully balance the column after previous metal removal, then use 0.1-0.3mol/L metal salt solution (chloride or sulfate soluble metal salts are suitable) to pass through the column for 5-10 CV to chelate the fresh metal ions, and finally use 5-10 CV of distilled water to remove unbound metal ions.

### 3.6 Cleaning-In-Place (CIP)

1. Removal of proteins adsorbed by ion exchange: first use 2 mol/L NaCl solution to backwash 2 to 3 CV, then rinse with 3 to 5 CV with distilled water.

2. Removal of precipitated proteins, hydrophobic proteins and lipids: first use 1.0 mol/L NaOH to backwash the column (50 cm/h, contact time 1-2 h), contact for more than 12 hours to remove endotoxin. Then rinse the chromatography column with 10 CV of equilibration buffer, and finally rinse the chromatography column with 5 CV of distilled water.

3. Removal of hydrophobic proteins and lipids: first wash the chromatography column with 5-10 CV of 30% isopropanol in reverse phase (contact time 15-20min), then rinse the chromatography with 10 CV of distilled water column.

#### 4. Storage

Sealed and stored at 4~30°C (preservation solution is 20% ethanol) in a ventilated, dry and clean place, do not freeze.

#### 5. Transportation

Avoid sunlight, rain, and heavy pressure during transportation. It is strictly forbidden to transport with toxic and hazardous materials.

#### 6. Precautions

6.1 The sample and chromatography media must be thoroughly equilibrated with equilibration buffer before column chromatography can be performed.

6.2 The loaded column bed must have a flat surface, with no channel flow or air bubbles, otherwise it should be reloaded.

6.3 During operation, ensure that the temperature of the column and the buffer are consistent to avoid air bubbles in the column bed, which affect the purification performance.

6.4 During the elution process, the flow rate should be strictly controlled.

6.5 During sample loading and the entire elution process, prevent the column surface from drying out.

#### 7. Ordering information

Product Name	References	Pack Size
Seplife® 6AG IDA/90	A4003012	25ml
	A4003013	100ml
	A4003014	500ml
	A4003015	1L
	A4003016	5L
	A4003017	10L

Production date: See label

Expiry date: 5 years, under proper storage conditions

**Manufacturer: Sunresin New Materials Co. Ltd.**

Add:No. 135, Jinye Rd, Xi'an Hi-tech Industrial Development Z

one, Shaanxi, 710076, China

[www.seplite.com](http://www.seplite.com)    [www.sunresin.com](http://www.sunresin.com)

E-mail: [info.lifescience@sunresin.com](mailto:info.lifescience@sunresin.com)

All information set forth herein is for informational purposes only. This information is general descriptive(introductory) information of SUNRESIN and its related products, technologies and services. Neither shall constitute the guarantee of SUNRESIN and its affiliates to products, technologies and services in specific fields and specific application conditions results, unless otherwise expressly noted. SUNRESIN and its affiliates assumes no obligation or liability for the information in this document. Customer is responsible for judging whether the information is appropriate for Customer's concrete demand and are obliged to understand whether the use of these products, technologies and services is permitted by the laws and regulations of their countries and relevant regions. Unless expressly stated, no freedom from infringement of use any patent or trademark or intellectual property rights owned by SUNRESIN or its affiliated companies under this document is to be inferred.